Lovibond® Water Testing

Tintometer® Group

Manual of Methods

MD 100 • MD 110 • MD 200

Chlorine

(EN) Manual of Methods

Page 4

(ES) Manual de Métodos

Página 16

(IT) Manuale dei Metodi

Pagina 28

(NL) Handboek Methoden

Zijde 40

(DE) Methodenhandbuch

Seite 10

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


Page 22

(PT) Métodos Manual

Página 34

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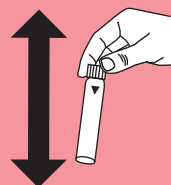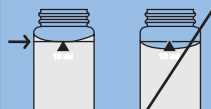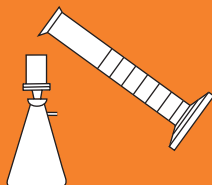Page 46

KS4.3 T / 20

Method name

Method number

Bar code for the detection of the methods

Measuring range

20

S:4.3

Display in the MD 100 / MD 110 / MD 200

Chemical Method

Instrument specific information

The test can be performed on the following devices. In addition, the required cuvette and the absorption range of the photometer are indicated.

Instrument Type	Cuvette	λ	Measuring Range
MD 200, MD 600, MD 610, MD 640, MultiDirect, PM 620, PM 630	ø 24 mm	610 nm	0.1 - 4 mmol/l $K_{S4.3}$
SpectroDirect, XD 7000, XD 7500	ø 24 mm	615 nm	0.1 - 4 mmol/l $K_{S4.3}$

Material

Required material (partly optional):

Reagents	Packaging Unit	Part Number
Alka-M-Photometer	Tablet / 100	513210BT
Alka-M-Photometer	Tablet / 250	513211BT

Application List

- Waste Water Treatment
- Drinking Water Treatment
- Raw Water Treatment

Notes

1. The terms Alkalinity-m, m-Value, total alkalinity and Acid demand to $K_{S4.3}$ are identical.
2. For accurate results, exactly 10 ml of water sample must be used for the test.

Language codes ISO 639-1

Revision status

EN Handbook of Methods 01/20

Performing test procedure

Implementation of the provision Acid capacity $K_{S_{4.3}}$ with Tablet

Select the method on the device

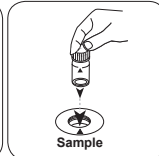
For this method, no ZERO measurements are to be carried out with the following devices: XD 7000, XD 7500

Fill 24 mm vial with **10 ml sample**.

Close vial(s).

Place **sample vial** in the sample chamber. • Pay attention to the positioning.

• • •

Dissolve tablet(s) by inverting.

Place **sample vial** in the sample chamber. • Pay attention to the positioning.

Press the **TEST (XD: START)** button.

The result in Acid Capacity $K_{S_{4.3}}$ appears on the display.

Chlorine HR (KI) T

M105

5 - 200 mg/L Cl₂

CLHr

KI / Acid

Material


EN

Required material (partly optional):


Reagents	Packaging Unit	Part Number
Chlorine HR (KI)	Tablet / 100	513000BT
Chlorine HR (KI)	Tablet / 250	513001BT
Acidifying GP	Tablet / 100	515480BT
Acidifying GP	Tablet / 250	515481BT
Set Chlorine HR (KI)/Acidifying GP 100 Pc. #	100 each	517721BT
Set Chlorine HR (KI)/Acidifying GP 250 Pc. #	250 each	517722BT
Chlorine HR (KI)	Tablet / 100	501210
Chlorine HR (KI)	Tablet / 250	501211

Determination of Chlorine HR (KI) with Tablet

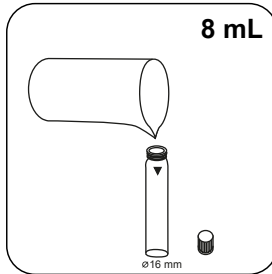
Select the method on the device.

Fill 16 mm vial with **8 mL sample**.

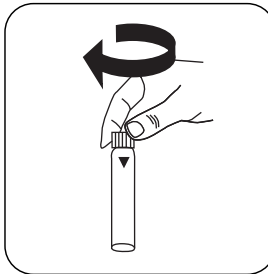
Close vial(s).

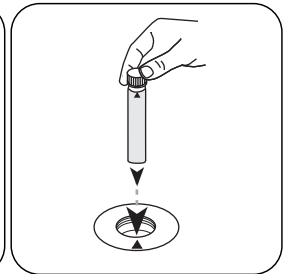Place **sample vial** in the sample chamber. • Pay attention to the positioning.

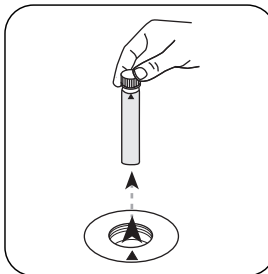
Press the **ZERO** button.


Remove **vial** from the sample chamber.

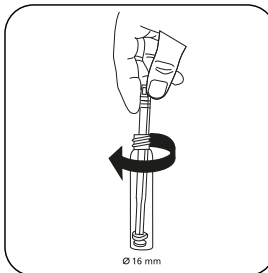
Add **Chlorine HR (KI) tablet**.


Crush tablet(s) by rotating slightly.

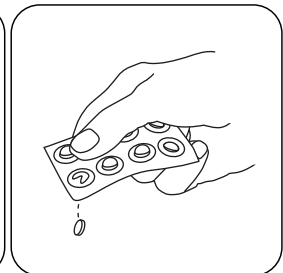
Add **ACIDIFYING GP tablet**.

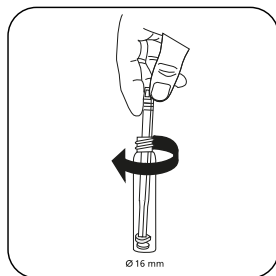
Crush tablet(s) by rotating slightly.

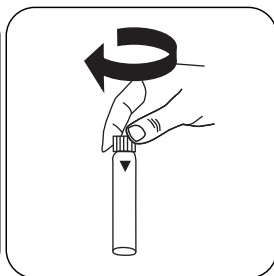
Close vial(s).

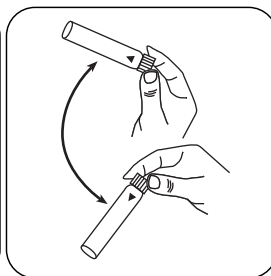Dissolve tablet(s) by inverting.

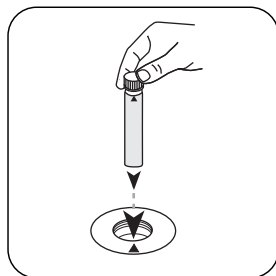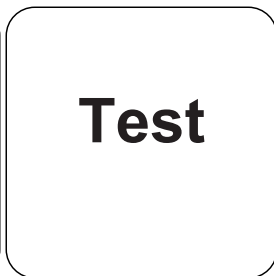Place **sample vial** in the sample chamber. • Pay attention to the positioning.

Press the **TEST** (XD: **START**) button.

The result in mg/L Chlorine appears on the display.

Chemical Method

KI / Acid

Appendix

Interferences

Persistent Interferences

- All oxidising agents in the samples react like chlorine, which leads to higher results.

Method Validation

Limit of Detection	1.29 mg/L
Limit of Quantification	3.86 mg/L
End of Measuring Range	200 mg/L
Sensitivity	83.96 mg/L / Abs
Confidence Intervall	1.14 mg/L
Standard Deviation	0.45 mg/L
Variation Coefficient	0.45 %

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Lovibond® and Tintometer® are Trademarks of
Tintometer Inc.