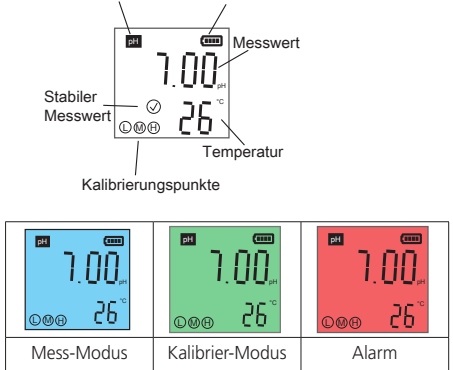
Display

Mess-Modus Kalibrier-Modus Alarm

Tastaturfunktion

Kurzer Tastendruck (1 Sek.) - Langer Tastendruck (3 Sek.)

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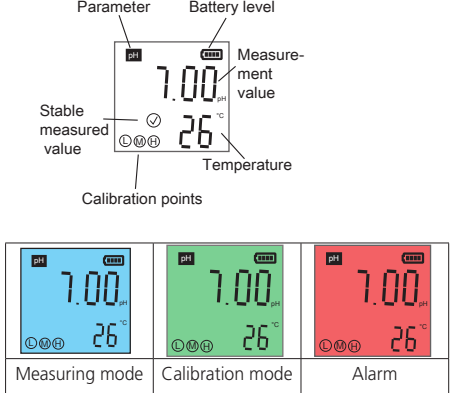
Display

Measuring mode Calibration mode Alarm

Keyboard function

Short keystroke (1 sec.) - Long keystroke (3 sec.)

Table with 3 columns: Action (Short/Long press), Mode (Measuring/Calibration), and Function (e.g., power on/off, backlight, settings).

Settings

In off mode, press 3 seconds: Opening Settings.


Navigate Confirm

Blue marked = factory setting

Table with 3 columns: Parameter (pH calibration series, alarm limit), Selection (USA, NIST, pH 0-14), and Meaning (description of the setting).

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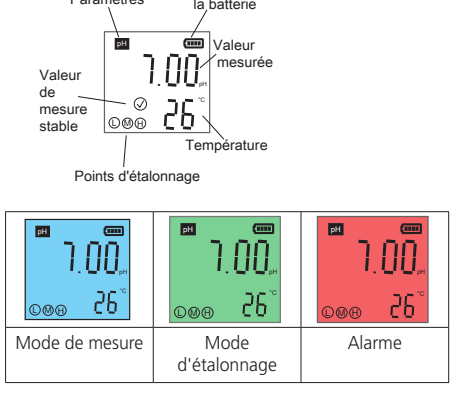pH Kalibrierung


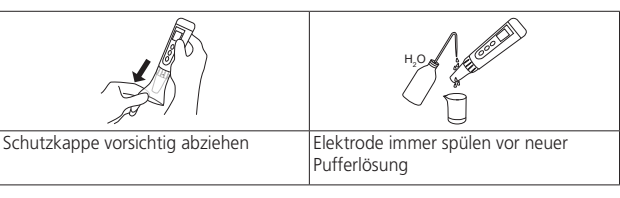
Table with 2 columns: 1-Punkt-Kalibrierung and 2-Punkt-Kalibrierung. Includes numbered steps and illustrations for each.

Table with 3 columns: Parameter (Auto-Hold, Backlight, temperature unit, factory calibration), Action (On/Off/1 min/Off/Yes/No), and Meaning (description of the setting).

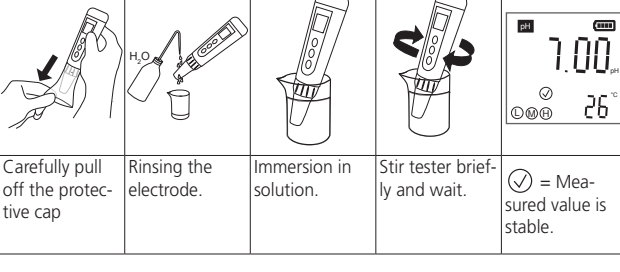pH measurement

Note: Press at "Hold" for new measurement recording.

pH Calibration


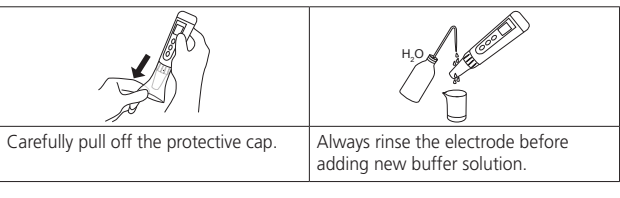
Table with 2 columns: 1-point calibration and 2-point calibration. Includes numbered steps and illustrations for each.

After successful 1, 2 or 3 point calibration, the display shows:

Table showing the display icons for 1-point, 2-point, and 3-point calibration.

Table with 3 columns: Parameter (Max. limit alarm, Auto-Hold, backlight, temperature unit, factory calibration), Action (On/Off/1 min/Off/Yes/No), and Meaning (description of the setting).

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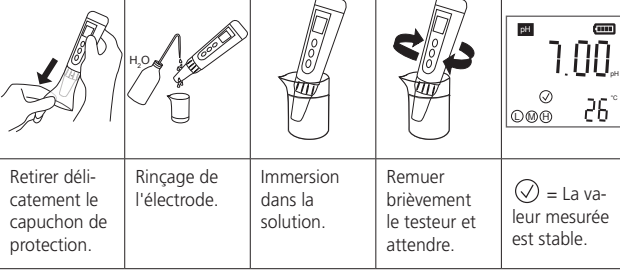
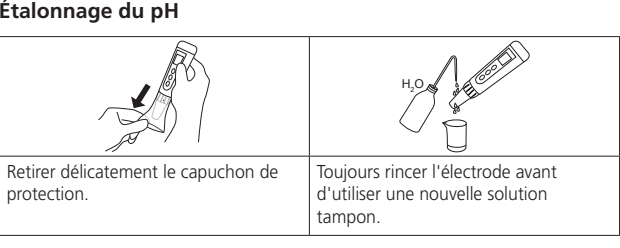
Table with 2 columns: Calibrage en 1 point and Calibrage à 2 points. Includes numbered steps and illustrations for each.

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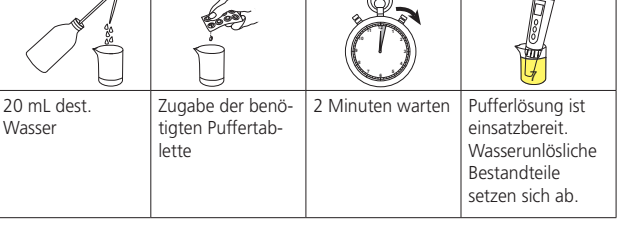Table with 2 columns: Parameter (pH, temperature, backlight, hold, auto-off, selection unit, housing protection, batteries) and Specification (range, resolution, accuracy, etc.).

Table with 2 columns: Austausch der Elektrode and Aufbewahrung der Elektrode. Includes illustrations for electrode replacement and storage.

Austausch Batterien


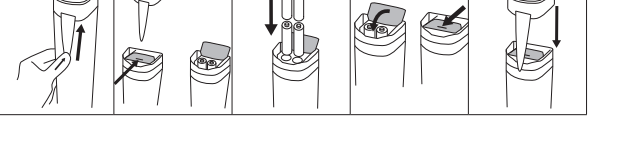
Table showing the display icons for 3-point calibration and pH 4/7/10.

Note: Other pH calibration series possible according to settings.

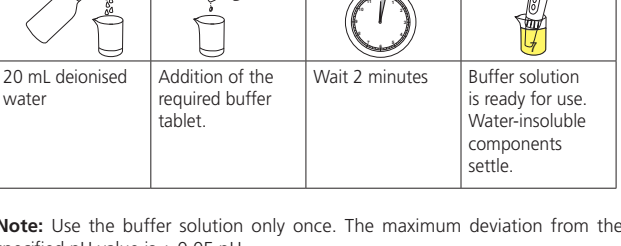Preparation of the buffer solution


Note: Use the buffer solution only once. The maximum deviation from the specified pH value is ± 0.05 pH.

SD 10 pH/T Specifications

Table with 2 columns: Parameter (pH, temperature, background light, hold, auto-off, selection unit, housing protection, batteries) and Specification (range, resolution, accuracy, etc.).

Table with 2 columns: Replacing the electrode and Storage of the electrode. Includes illustrations for electrode replacement and storage.

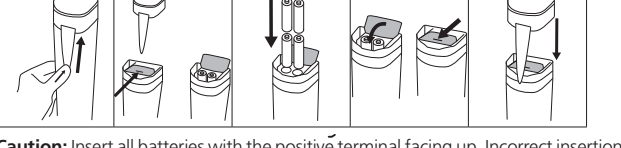Exchange batteries

Caution: Insert all batteries with the positive terminal facing up. Incorrect insertion of the batteries could result in damage to the meter.

Note: Make sure that the battery cover is completely closed to ensure water tightness.

Error messages and procedure

Table with 3 columns: Reason and Procedure for error messages.

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Zubehör

Table with 2 columns: Art.-Nr. (Article number) and Beschreibung (Description) of accessories.

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Table with 3 columns: Calibration mode, Error code (Err 1, Err 2), Reason, and Procedure.

Note: In case of heavy contamination, carefully remove the sensor protection and rinse the electrode.

Accessories

Table with 2 columns: Art.-No. (Article number) and Description of accessories.

Warranty

The warranty period for the SD10 pH is 2 years from the date of purchase and 6 months for the electrode.

Complete instructions with accessories list and further error messages can be found at www.lovbond.com (Download Centre).

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