CHECKIT[®] Comparator


pH (Bromthymolblau) 6.0 - 7.6

pH (Bromothymol Blue) 6.0 - 7.6

pH (Bleu de bromothymol) 6.0 - 7.6

pH (Azul de bromotimol) 6.0 - 7.6

pH (Blu di bromotimolo) 6.0 - 7.6

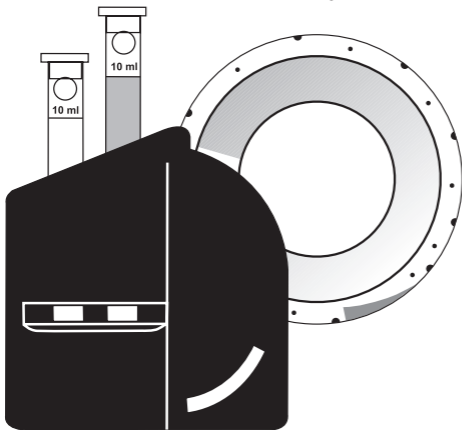DE

EN

FR

ES

IT

Hinweis:

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




Note:

Read the safety data sheets for the chemicals used. Use the recommended personal protective equipment and follow the disposal instructions.

Place the CHECKIT® Disc in the comparator with numeric values facing the operator.

**pH (Bromothymol Blue)**

1. Rinse the cuvettes with the water sample. Then fill both cuvettes with the water sample to the 10 mL mark. Close a cuvette with the lid and place it in the left chamber of the comparator as a blank. The dot on the cuvette must point to the viewer.
2. Add one BROMOTHYMOL BLUE tablet in the other cuvette. Do not touch the tablet with your fingers.
3. Close the cuvette with the lid and swirl slightly to dissolve the tablet faster.
4. After color development, place the cuvette in the right chamber of the comparator. The dot on the cuvette must point to the viewer.
5. Match the colour of the water sample with the CHECKIT® DISC.
The matching colour indicates the pH-value.

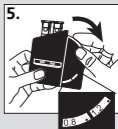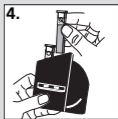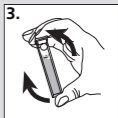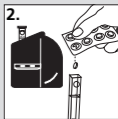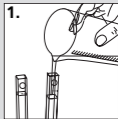Note:

pH values below 6.0 will always result in a yellow coloration.
pH values above 7.6 will always result in a blue coloration.
Water samples with low carbonate hardness (SBV $4.3 < 0.7$ mmol/l) may produce inaccurate pH values.

Important:

After each measurement, rinse the cuvettes thoroughly.
Store CHECKIT® Disc in the dark.
If stored properly, the color discs have a shelf life of 3 years from the date of receipt.

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